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09.06.11

AC 5125.25

This GM tube is suitable for measuring low levels of radiation where the small number of counts would give rise to large uncertainties. The GM tube has quite a large window, and will typically give 3 times higher count rates than a standard GM tube – for instance in half-life measurements with the isotope generator (no. 5112.00). This tube is also very well suited for detecting radon daughters.

As the tube's window is quite large – diameter 28.6 mm – it is very fragile. To minimize the risk of destroying the window, the tube is provided with a protective grating. This grating reduces the number of counts with approximately 15 %, but still the tube is superior to standard GM tubes. The grating is held by a threaded flange and may **with caution** be removed to obtain the maximum number of counts. (The window should face downwards during this operation.)

As extra protection the tube is provided with a protective cap. It is recommended always to use this cap for protection when the tube is not in use. Do not cover the small hole in the cap when it is mounted or removed.

The GM tube is provided with a BNC-plug.

Please notice that the GM-tube has no built-in anode resistor. Our GM counters are provided with a built-in anode resistor.

Technical specifications:

The tube is sensitive to alpha, beta and gamma radiation.

The length of the tube is 41 mm.

The window is retracted 7.5 mm from the front of the housing (without cap).

Plateau threshold voltage (V_{b1}): 450 V

Plateau length ($V_{b2} - V_{b1}$): 250 V

Recommended supply voltage: 500 V

Window specifications:

Thickness: 2.0 til 3.0 mg/cm²

Usefull diameter: 28.6 mm

Material: Mica

Gas filling: Neon and Halogen

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© A/S Søren Frederiksen, Ølgod