# Scout<sup>®</sup> STX Series Balances

## User Guide

EN

ES

FR

DE

IT

PT

HU

SV

DA

PL

CS

NL

# 1. SAFETY INFORMATION

## Definition of Signal Warnings and Symbols

- WARNING** For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
- CAUTION** For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
- Attention** For important information about the product.
- Note** For useful information about the product.

## Warning Symbols

General hazard

Electrical shock

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- Do not operate the equipment in hazardous or explosive environments.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

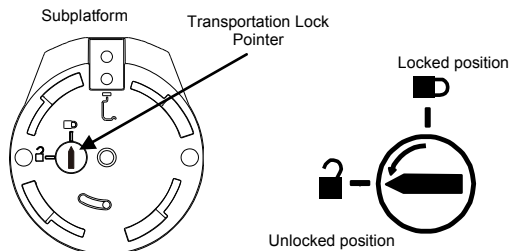
# 2. INSTALLATION

## 2.1 Selecting the location

The location must be sturdy, flat and level. Avoid locations with excessive air current, vibrations, heat sources or rapid temperature changes. Allow sufficient space around the instrument.


## 2.2 Releasing the Transportation Lock

On top of the balance, turn the pointer 90 degrees counter-clockwise to unlock.

### 2.3 Installing the Weighing Pan

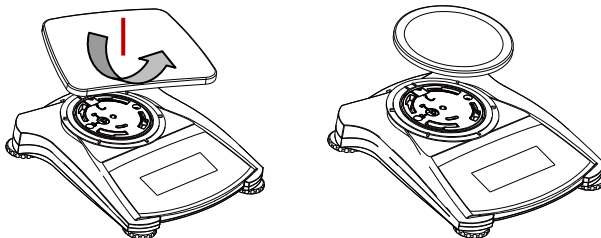Balances with a rectangular pan are placed into the sub-pan as shown and rotated counter-clockwise until it locks. Round pans are placed straight down on sub-pan.

### 2.4 Levelling the Instrument

To level the instrument, adjust the feet so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.

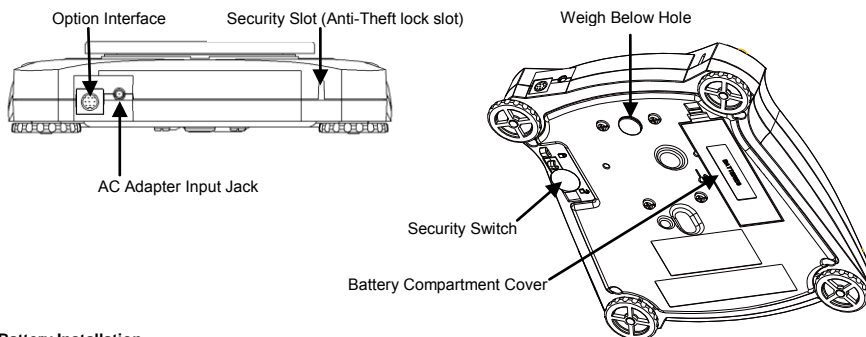
### 2.5 Connecting Power

#### AC Adapter Installation

Connect the AC adapter power cord to the instrument's power input connector, and then connect the AC plug to a suitable electrical outlet.

**Attention:** Only use a AC adapter specified by OHAUS.

#### Battery Installation

Install the four "AA" batteries with polarity as shown in the battery compartment.

### 2.6 Initial Calibration

For best results, the instrument's calibration should be checked prior to first use. If adjustment is needed, refer to the Calibration section of the instruction manual.

### 3. OPERATION

#### 3.1 Controls

| Button                          | ZERO  | TARE                                     |
|---------------------------------|---|--|
| Primary Function (Short Press)  | <b>On</b><br>Turns the balance on<br><b>Zero</b><br>If balance is On, sets Zero | <b>Tare</b><br>Enter/clear a tare value. |
| Secondary Function (Long Press) | <b>Off</b><br>Turns the balance Off.  |  |

#### 3.2 Overview of Display, Home Screen

This equipment utilizes a touch-sensitive display. *Touch* areas and Buttons to control the equipment's functions.

##### Main Application Screen

|   |  |   |
|---|--|---|
| <p>Application</p> <p>Instructional Messages</p> <p>Stability (*), Net (&gt;0&lt;) indicators</p> <p>Reference Fields</p> | <p>The screenshot shows the 'Dynamic Weighing' application. At the top, it displays 'Max 220 g d = 0.001 g' and the date/time '9/28/2015 11:33'. Below this is the instruction 'Place sample on the pan to begin averaging.' and a large result field showing '0.000 g'. At the bottom, there are three buttons: 'Item Settings', 'Print', and 'Menu'. Reference fields show 'Gross: 0.000 g', 'Averaging Time: 5 s', 'Tare: 0.000 g', and 'Automatic Mode: On'.</p> | <p>Capacity and readability</p> <p>Result Field:<br/>Information varies by application<br/>Touch <b>g</b> to change unit</p> <p>Application Buttons:<br/>Function vary by application</p> |
|---|--|---|

#### 3.3 MENU & SCREEN NAVIGATION

Touch **Menu** to open the menu list.

- 

**Calibration:**  
Touch to view calibration options.
- 

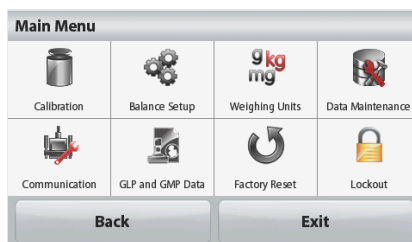
**Balance Setup:**  
Touch to view balance settings.
- 

**Weighing Units:**  
Touch to view weighing units.
- 

**Data Maintenance:**  
Touch to view data maintenance options.
- 

**Communication:**  
Touch to view COM Device Settings and Print Settings.
- 

**GLP and GMP Data:**  
Insert user data for traceability.

- 

**Factory Reset:**  
Touch to do a Factory reset.
- 

**Lockout:**  
Touch to access the Lockout menu list.

### 3.3 Using the Balance

#### Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Add sample to the pan or container. The display shows the weight of the sample.

## 4. MAINTENANCE

### 4.1 Calibration

When the Balance is first installed, and when it is moved to another location, it must be calibrated to ensure accurate weighing results.

### 4.2 Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

### 4.3 Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you.

## 5. TECHNICAL DATA

The technical data is valid under the following conditions:

- Indoor use only
- Altitude: Up to 2000 m
- Operating temperature range: +5 °C to +40 °C
- Specified temperature range: +10 °C to +40 °C
- Relative humidity: 10% to 80% at 31°C, decreasing linearly to 50% at 40°C, non-condensing
- Mains supply voltage fluctuations: up to ±10% of the nominal voltage
- Installation category II
- Pollution degree: 2

#### SPECIFICATIONS:

| MODEL                     | STX123   | STX223  | STX222 | STX422 | STX622 | STX1202 |
|---------------------------|--|---------|--------|--------|--------|---------|
| Capacity                  | 120 g  | 220 g   | 220 g  | 420 g  | 620 g  | 1000 g  |
| Readability d             | 0.001 g  | 0.001 g | 0.01 g | 0.01 g | 0.01 g | 0.01 g  |
| Repeatability (std. dev.) | 2 d  | 2 d     | 1 d    | 1 d    | 1 d    | 2 d     |
| Linearity                 | 3 d  | 3 d     | 1 d    | 1 d    | 2 d    | 3 d     |
| Power                     | AC power adaptor input 100-240 V 50/60 Hz and output 5V/1A DC, or 4 AA batteries |         |        |        |        |         |

| MODEL                     | STX2202   | STX421 | STX621 | STX2201 | STX6201 | STX8200 |
|---------------------------|---|--------|--------|---------|---------|---------|
| Capacity                  | 2200 g  | 420 g  | 620 g  | 2200 g  | 6200 g  | 8200 g  |
| Readability d             | 0.01 g  | 0.1 g  | 0.1 g  | 0.1 g   | 0.1 g   | 1 g     |
| Repeatability (std. dev.) | 2 d   | 1 d    | 1 d    | 1 d     | 1 d     | 1 d     |
| Linearity (g)             | 3 d   | 1 d    | 1 d    | 1 d     | 2 d     | 1 d     |
| Power                     | AC power adaptor input 100-240 V 50/60 Hz and output 5 V/1A DC, or 4 AA batteries |        |        |         |         |         |

## 6. COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

| Mark | Standard   |
|------|--|
|      | This product complies with the EU Directives 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .  |
|      | This product complies with the EU Directive 2002/96/EC (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> . |
|      | AS/NZS 61000.6.1, AS/NZS 61000.6.3   |

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P/N 30269065 B © 2015 Ohaus Corporation, All rights reserved
Ohaus Corporation